
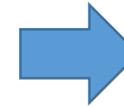
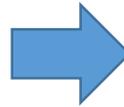


UAB Spill Kit Procedures

Absorbent Pad Spill Kit

Step One: Locate and don the appropriate PPE for your spill

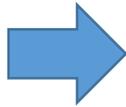


Step Two: Locate your labs Absorbent Spill Kit


Step Three: Take the absorbent barrier out of the kit and place around the spill.

Step Four: Place the absorbent pad directly on the spill. It may take more than one.

Step Five: Get the Hazardous Materials bag out of your spill kit. Place the used absorbent pads in the bag.

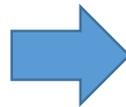

Step Six: Put an UAB Hazardous Material Label on the outside of the bag and properly fill it out.

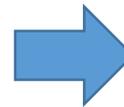


Step Seven: Make sure the bag is securely tied and ready for waste removal. Place in your lab's secured Satellite Accumulation Area.





UAB Spill Kit Procedures

Dry Spill Kit

Step One: Locate and don the appropriate PPE for your spill.

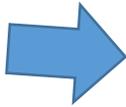

Step Two: Locate your labs Dry Spill Kit.

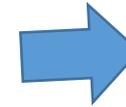
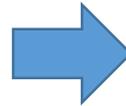


Step Four: Cover the spill completely with kitty litter



Step Six: Place the wet kitty litter in the appropriate waste container

Step Five: Once the spill has been absorbed, you need to sweep it up

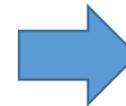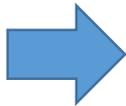

Step Seven: Make sure your waste container is properly labeled and ready for waste removal. Place in your labs Satellite Accumulation Area

